AGITADOR VORTEX RSLAB-6 / RSLAB-6PRO
VORTEX STIRRER RSLAB-6 / RSLAB-6PRO
AGITATEUR VORTEX RSLAB-6 / RSLAB-6PRO

Tel. +34 902 11 79 29
www.infoagro.com/tienda

RS Lab

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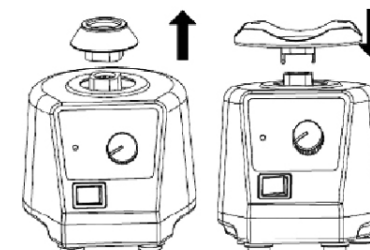
Spanish 2-7
 English..... 8-13
 French..... 14-19

INDEX OF CONTAINS

Preface and Warranty..... 8
 1. Safety instructions 9
 2. Proper use 9
 3. Inspection 9
 4. Trial run10
 5. Operation
 a. Continuous operating11
 b. Touch operating11
 6. Maintenance and cleaning11
 7. Associated standards and regulations12
 8. Technical data12
 9. Accessories13

Welcome to the vortex stirrer RSLAB-6 (without speed control) and RSLAB-6PRO (with speed control) user manual. Users should read this manual carefully, follow the instructions and procedures, and beware of all the cautions when using this instrument.

When help needed, you can always contact the service department of manufacturer or your supplier for technical support. Please provide the customer care representative with the following information: Serial number, description of problem, methods and procedures adopted to resolve the problems and your contact information.

These RSLAB vortex stirrers are warranted to be free from defects in materials and workmanship under normal use and service for a period of 12 months from the date of invoice. The warranty is extended only to the original purchaser. It shall not apply to any product or parts which have been damaged on account of improper installation, improper connections, misuse, accident or abnormal conditions of operation.

For claims under the warranty please contact your local dealer. You may also send the instrument direct to our works, enclosing the invoice copy and giving reasons for the claim.

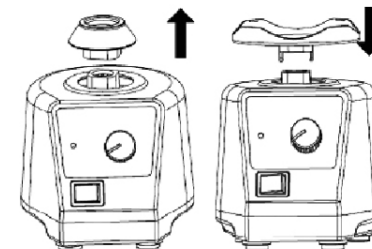
9- ACCESSORIES

- Code:** 90681581
 Standard top. For test tubes and small vessels up to Ø 30 mm
 Continuous operation / Touch operation
- Code :** 90681511
 Universal top 100 mm plate attachment with rubber pad
 Continuous operation
- Code :** 90681521
 Tube adapter, with 90681511. For 48 Eppen. tubes 6 mm
 Continuous operation
- Code :** 90681531
 Tube adapter, with 90681511. For 15 test tubes Ø 10 mm
 Continuous operation
- Code :** 90681541
 Tube adapter, with 90681511. For 12 test tubes Ø 12 mm
 Continuous operation
- Code :** 90681551
 Tube adapter, with 90681511. For 8 test tubes Ø 16 mm
 Continuous operation
- Code :** 90681561
 Tube adapter, with 90681511. For 8 test tubes Ø 20 mm
 Continuous operation
- Code :** 90681591
 Platform pad, with 90681511. For < Ø 99mm tubes and small vessel
 Touch operation
- Code** 90681582
 Spare rubber foot For RSLAB-6PRO

Note: 90681511 at 90681561 accessories are only suitable with the adjustable speed model (RSLAB-6PRO)

Serial number and pattern number and accessories names on the type plate must be supplied when ordering other accessories.

Changing attachment:

The instrument is a sophisticated electronic product, avoiding severe shake in work and transportation. The instrument must be cleaned and put in into the initial packaging carton before sending to service for repair, avoiding the contamination of hazardous.

Switch off the instrument and put it in the dry, clean, stable place at room temperature in long-term disuse.

7- ASSOCIATED STANDARDS AND REGULATIONS

Construction in accordance with the following safety standards:
EN61010-1

Construction in accordance with the following EMC standards:
EN61326-1

Associated EU guidelines:

EMC-guidelines: 89/336/EWG

Instrument guidelines: 73/023/EWG

8- TECHNICAL DATA

<i>Voltage</i>	220-230/110-120 VAC
<i>Frequency</i>	50/60 Hz
<i>Power</i>	60 W
<i>Shaking movement</i>	Orbital
<i>Orbital diameter</i>	4 mm
<i>Motor type input</i>	58 W
<i>Motor type output</i>	10 W
<i>Permissible ON time</i>	100 %
<i>Speed range</i>	0-2500 rpm (adjustable) 2500 rpm (fixed)
<i>Speed display</i>	Scale
<i>Run type</i>	Continuous / Touch operation
<i>Dimensions</i>	127x130x160 mm
<i>Weight</i>	3,5 kg
<i>Permitted ambient temperature</i>	5-40 °C
<i>Permitted relative humidity</i>	80 %
<i>Protection class acc. to DIN 60529</i>	IP 21

1- SAFETY INSTRUCTIONS

Warning!

*Read the operating instructions carefully before use the instrument.
Ensure that only trained staff work with the instrument*

Protective ground contact!

- ◆ Make sure that socket is earthed (protective ground contact) before use.
- ◆ When work, wear the personal guard to avoid the risk from splashing and evaporation of liquids and release of toxic or combustible gases.
- ◆ Keep medium vessels vertical as much as possible, mix the medium in the minimum pressure required and use attachment supplied when mixing a mass of different medium.
- ◆ Set up the instrument in a spacious area on a stable, clean, non-slip, dry and fireproof surface, do not operate the instrument in explosive atmospheres, with hazardous substances or under water.
- ◆ Gradually increase the speed, reduce the speed if the stirring bar breakaway because of too high speed and if the instrument is not running smoothly, or if the container moves on the stage
- ◆ Check the instrument and accessories before handle it for damage each time you use them. Do no use damaged components. Safe operation is only guaranteed with the accessories described in "Accessories" chapter. Accessories must be securely attached to the instrument and cannot come off by themselves. Always disconnect the plug before fitting accessories.
- ◆ Place a vessel at the centre of the mixing head and spread out others uniformly. A sharp edged vessel can wear out the mixing head.
- ◆ The instrument may heat up when in use.
- ◆ Protect the instruments and accessories from bumps and impacts.
- ◆ The voltage started on the label must correspond to the main power supply.
- ◆ The instrument may only be opened by qualified and trained technician

2- PROPER USE

The instrument is designed for mixing liquids in schools, laboratories or factories. This device is not suitable for using residential areas or other areas that may cause danger to the user or instrument as mentioned in chapter 1.

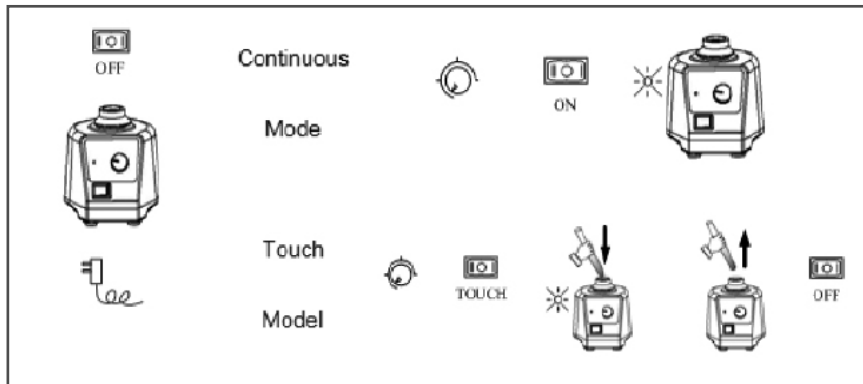
3- INSPECTION

- ◆ Unpack the equipment carefully and check for any damages which may have arisen during transport. If it happens, please contact manufacturer for technical support.

If there is any apparent damage to the system, please do not plug it into the power line

- ◆ The RSLAB-6 y RSLAB-6PRO includes the following items:
 - Main unit: 1 unit
 - User manual:1 unit
- ◆ Please retain all packing materials. Additional shipping charges will be assessed if packing is unavailable for the system warranty service.


4- TRIAL RUN

Ensure the two-direction switch must be in the middle (off state) before powering on.

If these operations above are normal, the instrument is ready to operate following the "Operation" chapter. If these operations are not normal, the instrument may be damaged. Please contact manufacturer for technical support.

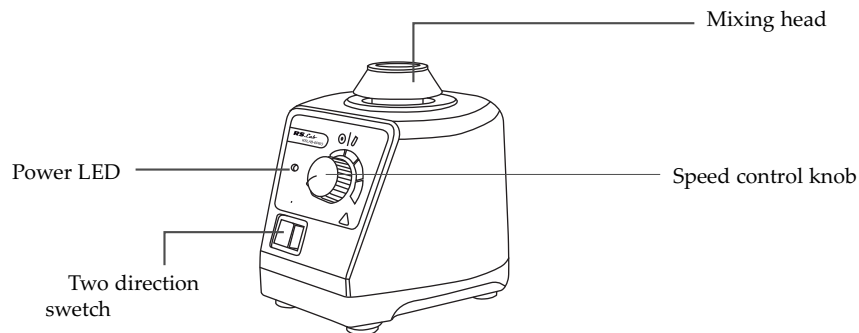5- OPERATION

1. Continuous operating

- ◆ Ensure the speed control knob to the lowest position. Put the instrument on the stable and safe place and plug in the main power.
- ◆ Push the two direction switch leftward to on and power LED is lit. The instrument begins to work.
- ◆ Turn the speed control knob to set the rated speed (adjustable model).
- ◆ If use the nonstandard accessories, please ensure the speed in indicate range (adjustable model).
- ◆ Push the two-direction switch rightward gently to off and LED power is closed and the instrument is turned off.

2. Touch operating

- ◆ Ensure the speed control knob to the lowest position. Put the instrument on the stable and safe place and plug in the main power.
- ◆ Push the two-direction switch rightward to the touch mode.
- ◆ Turn the speed control knob to set the rated speed (adjustable model).
- ◆ If a test tube is pressed into the mixing head vertically, the instrument will begin to work and the power LED will be lit.
- ◆ Push the two-direction switch leftward gently to off and the instrument is turned off.
- ◆ Place vessels such a test tube as vertical as possible into the centre of the mixing head in order to mixing smoothly.
- ◆ Adjust the speed control knob slowly in order to have the instrument work smoothly.

6- MAINTENANCE AND CLEANING

Proper maintenance can keep instruments working in a good state and lengthen its life-time.

Do not spray cleanser into the instrument when cleaning.

Unplug the power line when cleaning.

Only use cleanser that we advised as below:

- Dyes: isopropyl alcohol
- Construction materials: water containing tenside / Isopropyl alcohol
- Cosmetics: water containing tenside / Isopropyl alcohol
- Foodstuffs: water containing tenside
- Fuels: water containing tenside

Before using other method for cleaning or decontamination, the user must ascertain with the manufacturer that this method does no destroy the instrument.

Wear proper protective gloves during cleaning of the instrument

Electrical instruments may not be placed in the cleansing agent for the purpose of cleaning.